

## PerfeCta® NGS Quantification Kit for Illumina® Sequencing Platforms

Cat. Nos.	95154-500	PerfeCta NGS Quantification Kit – Illumina	Size:	500 x 20- $\mu$ L reactions	<b>Store at -25°C to -15°C protected from light</b>
	95155-500	PerfeCta NGS Quantification Kit – Illumina / ROX		500 x 20- $\mu$ L reactions	
	95156-500	PerfeCta NGS Quantification Kit – Illumina / Low ROX		500 x 20- $\mu$ L reactions	

### Description

Accurate quantification of the number of amplifiable library molecules is one of the most important factors for obtaining high quality read data with next-generation sequencing technologies. The PerfeCta NGS Library Quantification Kit uses real-time PCR to specifically quantify library molecules that possess the appropriate adapter tag at each end. These are the suitable template molecules for the bridge PCR used for Illumina NGS platforms. The included stabilized, pre-diluted DNA standards and pre-qualified primer set simplifies the library quantification process, minimizes pipetting errors, and ensures reproducible and precise qPCR results, even with dilute samples. The robust qPCR performance of PerfeCta SYBR® Green SuperMix provides accurate quantification of NGS libraries with varying fragment sizes or GC content.

### Instrument Compatibility

Different real-time PCR systems employ different strategies for the normalization of fluorescent signals and correction of well-to-well optical variations. It is critical to match the appropriate qPCR reagent to your specific instrument. Please consult the following table, or visit our web site at [www.quantabio.com](http://www.quantabio.com) to find the optimal kit for your instrument platform.

NGS Quant. Cat. No.	qPCR Reagent	Compatible Real-Time PCR Systems
95154-500	PerfeCta SYBR Green SuperMix	Bio-Rad CFX96™, CFX384™, Opticon™, MiniOpticon™, Chromo4™, iQ™5*, MyiQ™*, iCycler iQ®* (*requires persistent well factors) Cepheid Smart Cycler®, Qiagen/Corbett Rotor-Gene® Illumina® Eco™; Eppendorf Mastercycler® ep realplex Roche Applied Science LightCycler® 480, 96
95155-500	PerfeCta SYBR Green SuperMix, ROX	Applied Biosystems 7000, 7300, 7700, 7900, 7900HT, 7900HT Fast, StepOne™, StepOnePlus™
95156-500	PerfeCta SYBR Green SuperMix, Low ROX	Applied Biosystems 7500, 7500 Fast, ViiA™ 7, QuantStudio™ Agilent/Stratagene MX4000™, MX3005P™, MX3000P™

### Components

Name	Description	Part Number	Volume
Library Dilution Buffer (10x)	10X concentrated buffer containing 0.1M Tris-HCl (pH 8.0), 1 mM EDTA, and stabilizers	84287	2 x 3 mL
Illumina Primer Mix	Mix of 10 $\mu$ M (each) Illumina "P5" and "P7" primer sequences for Illumina sequencing libraries in 10 mM Tris-HCl (pH 8.0), 0.1 mM EDTA. Illumina forward primer: 5'-AAT GAT ACG GCG ACC ACC GA-3' Illumina reverse primer: 5'-CAA GCA GAA GAC GGC ATA CGA-3'	84280	0.3 mL
Illumina Standard 1	0.0005 pM linear, dsDNA standard in 1X library dilution buffer	84275	0.1 mL
Illumina Standard 2	0.005 pM linear, dsDNA standard in 1X library dilution buffer	84276	0.1 mL
Illumina Standard 3	0.05 pM linear, dsDNA standard in 1X library dilution buffer	84277	0.1 mL
Illumina Standard 4	0.5 pM linear, dsDNA standard in 1X library dilution buffer	84278	0.1 mL
Illumina Standard 5	5 pM linear, dsDNA standard in 1X library dilution buffer	84280	0.1 mL
qPCR Reagent	2X concentrated master mix for SYBR Green qPCR		
95154-500	PerfeCta SYBR Green SuperMix	2 x 95054-100	4 x 1.25 mL
95155-500	PerfeCta SYBR Green SuperMix ROX	2 x 95055-100	4 x 1.25 mL
95156-500	PerfeCta SYBR Green SuperMix Low ROX	2 x 95056-100	4 x 1.25 mL

### Storage and Stability

Store components in a constant temperature freezer at -25°C to -15°C protected from light upon receipt.

Repeated freezing and thawing will not affect product performance.

For lot specific expiry date, refer to package label, Certificate of Analysis or Product Specification Form

### Quality Control

Kit components are free of contaminating DNase and RNase. qPCR of the DNA Standards are assayed in triplicate reactions with PerfeCta SYBR Green SuperMix and must generate a standard curve with  $r^2 \geq 0.995$  and a PCR efficiency  $\geq 90\%$ . Dissociation curve analysis must present as a single melt peak with the expected  $T_m$ .

### NGS Library Sample Dilutions

Prepare 1X library dilution buffer with nuclease-free, molecular biology grade water. Do not use DEPC-treated water. Do not use autoclaved water. Properly maintained ultra-filtration systems generally deliver water of suitable quality.

Add 0.1 mL of 10X Library Dilution Buffer to 0.9 mL of nuclease-free water. Mix by vortexing. Larger or smaller quantities of 1X dilution buffer may be prepared depending on your experimental requirements. Each library sample requires approximately 0.3 mL. The 1X dilution buffer is stable at 2-8°C for 6 months.

The optimal dilutions for qPCR may vary depending on the nature of your NGS sequencing operations. Use the following example as a general guideline. It is important for one of your library dilutions to fall within the linear dynamic range of the supplied DNA standards.

- Prepare a 1:20 dilution of the library sample:  
Add 2 µL of library sample to a 1.5-mL tube containing 38 µL of 1X dilution buffer  
Mix by vortexing, and then centrifuge to collect contents
- Prepare 2 successive dilutions for qPCR as follows:  
1:2000 diln.                    Transfer 2 µL of the 1:20 dilution to a 1.5-mL tube containing 198 µL of 1X dilution buffer.  
    Mix by vortexing and then centrifuge to collect contents  
1:20,000 diln.                Transfer 5 µL of the 1:2000 dilution to a 1.5-mL tube containing 45 µL of 1X dilution buffer.  
Mix by vortexing and then centrifuge to collect contents

### qPCR Reaction Assembly

Preparation of a reaction cocktail is recommended to reduce pipetting errors and maximize assay precision. Use the appropriate PerfeCTa SYBR Green SuperMix for your qPCR instrument (see *Instrument Compatibility* section).

- Prepare the reaction cocktail:

Component	Volume for complete 20-µL rxn.	Volume for a 30-rxn reaction cocktail	Final Concentration
PerfeCTa SYBR Green SuperMix (2X)	10 µL	300 µL	1x
Illumina Primer Mix (10 µM each)	0.6 µL	18 µL	300 nM (each)
Nuclease-free water	5.4 µL	162 µL	
Template DNA	<u>4 µL</u>	_____	

**Note:** For smaller, or larger, reaction volumes scale all components proportionally. For example: when working with a 384-well plate carry out reactions in 10-µL (final) volumes using 8 µL of reaction cocktail and 2 µL of DNA standard or NGS library dilution.

Final Volume	20 µL	480 µL
--------------	-------	--------

Vortex gently to mix contents. Centrifuge briefly to collect components at the bottom of the reaction tube.

- Label 8 x 1.5-mL tubes A to H. Dispense 56 µL of reaction cocktail into each tube.
- Add 14 µL of the appropriate standard or NGS library dilution (1:2,000 or 1:20,000) to each tube. (See suggested plate loading below)
- Mix each tube by gently vortexing and centrifuge to collect contents.
- Dispense 20 µL from each tube into 3 wells for triplicate reactions for each DNA standard concentration or library dilution sample.

Row A: 3 x 20 µL of 1X dilution buffer	(NTC)
Row B: 3 x 20 µL tube "B": Standard 1	(0.0005 pM)
Row C: 3 x 20 µL tube "C": Standard 2	(0.005 pM)
Row D: 3 x 20 µL tube "D": Standard 3	(0.05 pM)
Row E: 3 x 20 µL tube "E": Standard 4	(0.5 pM)
Row F: 3 x 20 µL tube "F": Standard 5	(5 pM)
Row G: 3 x 20 µL tube "G": 1:2,000 Library Dilution	
Row H: 3 x 20 µL tube "H": 1:20,000 Library Dilution	
- Seal the plate and briefly centrifuge to collect contents to the bottom of each well

**Note:** Alternately, reaction cocktail (without DNA template) can be dispensed into the reaction plate (16 µL / well), and then DNA template added to each well to complete each qPCR reaction. Mix reactions after adding DNA template before PCR amplification.

## qPCR Amplification

1. Program your real-time qPCR instrument – reaction plate well assignments.
  - a. Define wells containing the DNA standards as “standards” with the appropriate concentration listed above.
  - b. Define NGS library dilution samples as “unknowns”
  - c. Group replicate reactions to allow the qPCR data analysis program to calculate average C<sub>q</sub> and concentrations based on the DNA standards.
2. Incubate the reaction plate in a real-time qPCR system as follows:

Initial denaturation:	95°C, 3 min
PCR cycling (35 cycles)	Denaturation: 95°C, 10 to 15 s
	Annealing: 60°C, 20s
	Extension: 72°C, 45s (collect and analyze data)
Dissociation (melt) curve	refer to your qPCR instrument manual ( <i>optional</i> )

Full activation of the hot-start DNA polymerase occurs within 30 seconds at 95°C. Initial denaturation times greater than 3 minutes are usually not required. However, amplification of GC-rich libraries may benefit from a longer initial denaturation step (5-10 min) to fully denature the template and minimize the potential for renaturation of highly stable and/or repetitive sequence regions. Increased extension time (1 min) may improve performance for libraries with longer (up to 1 kb) average insert sizes. For AT-rich libraries, use a 2-step cycling protocol with a combined annealing and extension step of 60°C for 1 min.

## Data Analysis

The reaction plate should be defined, or annotated, using the stated concentrations for each DNA standard. Since equal volumes of each DNA Standard and library dilution sample were used in each qPCR, a simple trend analysis of the standard curve (log pM concentration vs C<sub>q</sub> value) linear regression can be used to calculate the pM concentration for each library sample. This should be carried out automatically using the absolute (or standard curve) quantification algorithm provided with your qPCR detection system. There is no need to determine the template copy number or template concentration of each qPCR. Please consult your qPCR system documentation.

The Illumina DNA standards and primer mix generate a 426-bp amplicon (48.8% GC). A common problem with some NGS library quantification protocols is the use of DNA standards that are overly concentrated and generate qPCR data that are outside of the linear dynamic range for most qPCR instruments. Improper baseline settings result in compressions between the highest concentrated DNA standards, which in turn give rise to inflated PCR efficiencies and inaccurate library quantification results. The NGS DNA standards supplied with the PerfeCtа NGS Library Quantification Kits have been selected to avoid these artifacts and produce absolute quantification standard curves with exceptionally high linear regression correlation coefficients.

Size-adjusted concentration (pM) for each library dilution sample is obtained as follows:

$$\text{average concentration (pM)} \times (426 / \text{average fragment length})$$

The size-adjusted concentration of the original library sample is obtained by multiplying the value above by the dilution factor used for the qPCR (either 2,000 or 20,000). It is important to only use C<sub>q</sub> data from library dilution samples that fall within the linear dynamic range of the DNA Standards.

Use the size-adjusted concentration of the original library to determine the appropriate dilution to use for loading the flow cell for the bridge amplification. The optimal loading concentration may vary for different sample types or library construction methods, but in general, 5 to 10 pM generates an appropriate cluster density.

## Limited Label Licenses

Use of this product signifies the agreement of any purchaser or user of the product to the following terms:

1. The product may be used solely in accordance with the protocols provided with the product and this manual and for use with components contained in the kit only. QIAGEN Beverly, Inc. grants no license under any of its intellectual property to use or incorporate the enclosed components of this kit with any components not included within this kit except as described in the protocols provided with the product, this manual, and additional protocols available at [www.quantabio.com](http://www.quantabio.com). Some of these additional protocols have been provided by Quantabio product users. These protocols have not been thoroughly tested or optimized by QIAGEN Beverly, Inc.. QIAGEN Beverly, Inc. neither guarantees them nor warrants that they do not infringe the rights of third-parties.
2. Other than expressly stated licenses, QIAGEN Beverly, Inc. makes no warranty that this kit and/or its use(s) do not infringe the rights of third-parties.
3. This kit and its components are licensed for one-time use and may not be reused, refurbished, or resold.
4. QIAGEN Beverly, Inc. specifically disclaims any other licenses, expressed or implied other than those expressly stated.
5. The purchaser and user of the kit agree not to take or permit anyone else to take any steps that could lead to or facilitate any acts prohibited above. QIAGEN Beverly, Inc. may enforce the prohibitions of this Limited License Agreement in any Court, and shall recover all its investigative and Court costs, including attorney fees, in any action to enforce this Limited License Agreement or any of its intellectual property rights relating to the kit and/or its components.

©2018 QIAGEN Beverly Inc. 100 Cummings Center Suite 407J Beverly, MA 01915

Quantabio brand products are manufactured by QIAGEN, Beverly Inc.

Intended for molecular biology applications. This product is not intended for the diagnosis, prevention or treatment of a disease.



## Related Products

### sparQ DNA Library Prep Kit

Cat. No.	95191-024	Size:	24 reactions
	95191-096		96 reactions

### sparQ Adapter Barcodes Set A

Cat. No.	95193-A96	Size:	96 reactions
----------	-----------	-------	--------------

### sparQ Adapter Barcodes Set B

Cat. No.	95193-B96	Size:	96 reactions
----------	-----------	-------	--------------

### sparQ HiFi PCR Master Mix

Cat. No.	95192-050	Size:	50 reactions
	95192-250		250 reactions

The sparQ HiFi PCR Master Mix is a high efficiency, high-fidelity, and low bias PCR master mix for NGS workflows requiring DNA library amplification prior to sequencing. The included primer mix allows amplification of DNA libraries flanked by adapters containing the P5 and P7 Illumina® flow cell sequences.

## Trademarks

PerfeCTa is a registered trademark of QIAGEN Beverly Inc. SYBR is a registered trademark of Molecular Probes, Inc. Illumina is a registered trademark of Illumina, Inc. Applied Biosystems, StepOne, StepOnePlus, ViiA, QuantStudio and ROX are trademarks Life Technologies Corporation. Stratagene, MX3000P, MX3005P and MX4000 are trademarks of Agilent Technologies, Inc. LightCycler is a registered Trademark of Roche. Mastercycler is a trademark of Eppendorf. Rotor-Gene is a registered trademark of Qiagen GmbH. SmartCycler is a trademark of Cepheid. CFX96, CFX384, iCycler iQ, iQ5, MyiQ, Opticon, MiniOpticon and Chromo4 are trademarks of Bio-Rad Laboratories.