

## Plugging in Q qPCR Cycler

Please follow the steps below to ensure proper use of your Q qPCR Cycler.

**MAKE SURE POWER PACK IS NOT CONNECTED TO MAINS POWER WHEN CONNECTING TO Q qPCR Cycler**

Plug the power cord into the adaptor and insert the adaptor into the back of the instrument.

---

**CAUTION**



To avoid electrical arcing, ensure that the power adapter is not plugged into a wall socket before plugging into the instrument.

---

**Plug the power cord into a wall socket and switch the power on at the socket.**

**Power the instrument 'On' using the power switch at the back of the instrument.**

An illuminated blue light at the front of the instrument will show the instrument is powered on.